# 동서과학(주) Compact Pump Drive MANUAL

TC 600

## CONTENTS

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

## 11. COMM SETTING ( RS485 communication setting )

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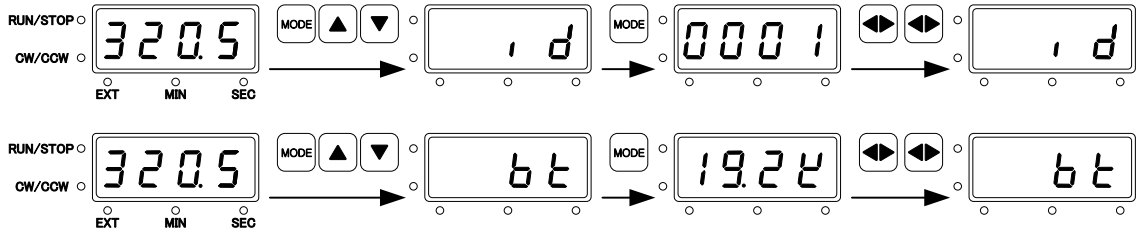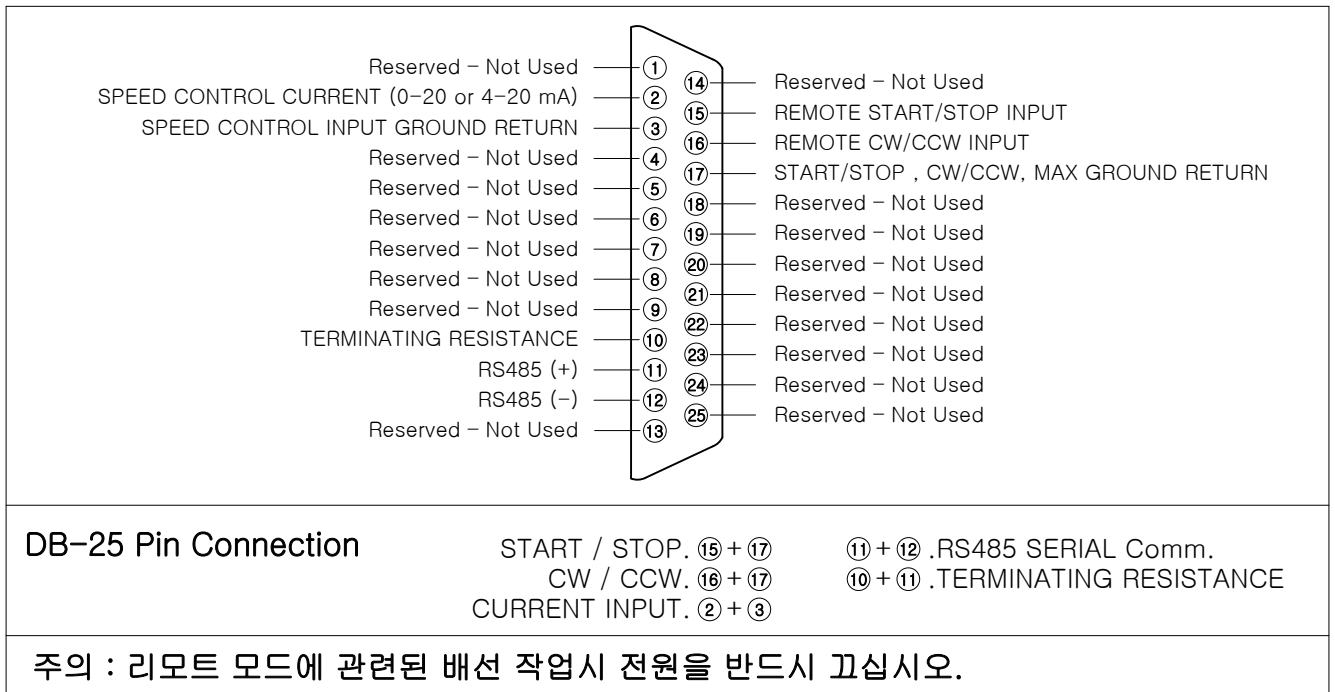## 14. SPECIFICATION

Description		MODEL	
Product		Compact Pump Drive	
Output		TC 600	
	Speed	0.1 – 600 rpm/min	
	Torque Output	1.5 ~ 30 (kgf – cm)	
	Speed accuracy	±0.1%	
	Display	7 Segment LED	
Input			
	Supply voltage	230V~, 50/60Hz	
	Fuse	T3.0A 250VAC	
	Current	1.0A-230Vrms	
Remote Input			
		START/STOP, CW/CCW (contact closure )	
	Current input	04-20mA,250Ω, ±0.5% full scale	
Construction			
	Dimension	180W x 270L x 221H (mm)	
	Weight	6.6 kg	
	Enclosure	ABS V0, Aluminum paint coated	
Environment			
	Temperature, Operating	0° to 40°C (32° to 104°F)	
	Temperature, Storage	-25° to 65°C (-13° to 149°F)	
	Humidity (non-condensing)	10% to 90%	
	Altitude	Less than 2000m	
	Pollution Degree	Pollution Degree 2 (Indoor use-lab, office)	
		MAINS supply voltage fluctuations up to ±10 % of the nominal voltage	
		Ptransient overvoltages typically present on the MAINS supply. “Installation Category II” or Value of transient voltage.	

## 13. CAUTION

**DANGER** : High pressure is flowing and can cause an electric shock.  
Be careful when you attempt to repair the inside of the machine.

**WARNINGS** : Tubing breakage may cause fluid to spray.  
Protect the driver and equipment by taking appropriate action on the pump drive.

To prevent your fingers or loose clothing from getting caught in the drive system, turn off the drive before removing or installing the tubing.

**CAUTION** : When connecting the external remote control cable to the pump drive, turn off the power to avoid damaging the pump drive.

**CAUTION** : Use the supplied authentic products for pump drive, pumphead and tubing. Failure to comply with the above may cause the tubing to break and leak when using the pump drive.

**CAUTION** : To avoid electric shock, make sure to ground the machine using a grounded power cord. Do not operate in humid places.

**CAUTION** : To prevent your fingers or loose clothing from getting caught in the drive system, turn off the drive before removing or installing the tubing.

**CAUTION** : Refer to the user manual to reduce the risks and ensure correct device operation without dangerous situations.

**CAUTION** : The surface is hot, so don't touch it.

**This product is not designed for medical using.  
Refer to the manual for user safety and device safety.**

14. The warranty lasts one year and excludes careless use or expendable parts.

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